



**Pneumatic Scale**  
**Angelus**

**CARR Centritech**  
**Separation Systems**

## Powerfuge® P12



### Features

- Drier solids and clear liquids
- Sealed automatic operation
- Full CIP/SIP designs
- cGMP construction
- Designs available for ASME-rated boundary, hazardous locations and aseptic operations
- State-of-the-art manufacturing facility located in Clearwater, Florida USA

## Powerfuge® P12

PneumaticScaleAngelus' unique Powerfuge® P12 System provides continuous flow two-phase separation of solids and liquids at rates up to 500 liters per hour for batches up to 4,000 liters. Using an innovative design, the feed is introduced through a stationary pipe and accelerated to full rotational speed before entering the titanium-alloy bowl. In this system, centrifugal forces as 20,000g permit separation of even sub-micron particles. Compressed solids are periodically removed with a fully automated scraping cycle. After the solids have been discharged, automated CIP/SIP cycles can be performed via the PLC controlled operating sequence.

The Powerfuge® P12 is designed for ease operation, with seal feed and drain lines allowing for aerosol-free operation. Similarly, no special tools are required to disassemble the system. Integrated turnkey systems incorporating feed delivery options, CIP and SIP skids, temperature control, and closed-loop centrate pumping under vacuum are available.

The Powerfuge® P12 system meets cGMP and CE requirements and has been validated to meet chemical and pharmaceutical industry protocols. The P12 is engineered to optimize ease of operation while maximizing throughput; each step in the process can be readily automated so that the time spent on routine maintenance can be minimized.

For more information, please contact us at:

**CellD**  
20bis, rue du Chapitre  
F-30150 ROQUEMAURE  
Tel : +33 (0)4 66 82 82 60  
Fax : +33 (0)4 66 90 21 10  
Email : [contact-info@celld.com](mailto:contact-info@celld.com) [www.celld.com](http://www.celld.com)

